

SOP 051 - Sequence Setup - Thermo Xcalibur

- 5.1.27.2. In the next screen, click directly on the “Yes” under the “Start Instrument” column next to the incorrect instrument.
- 5.1.27.3. Click again under the “Start Instrument” column next to the correct instrument so that “Yes” is displayed (ask the Instrument Chemist for additional help if needed)
- 5.1.27.4. Click “OK” and go to step 5.1.27.
- 5.1.28. Also ensure that the “Run Rows” field contains the correct starting and ending rows for the sequence. If not, enter the correct rows.
- 5.1.29. Click OK to begin the run.
- 5.1.30. Stay by the instrument to ensure that the first sample is injected properly before leaving.

6. References

- 6.1. Thermo Scientific. *Thermo Xcalibur Acquisition and Processing User Guide*. Sept. 2010. User Manual. USA.

Uncontrolled Copy