CHARLOTTE-MECKLENBURG POLICE DEPARTMENT

Crime Laboratory – Biology Section Biology QAM 8



Certificate of Quality Control Automate Performance Check

Specifications:

The Automate will be checked annually and after any repair or service with a known and a blank sample. Upon QPCR, the known must contain at least 0.075ng/µl of DNA and the blank must not have any DNA. If DNA is detected in the QPCR results for the blank, the test must be repeated by a DNA Team Leader or the Chief Criminalist.

Reason for the performance check:	Ann	ual
After repair, service, or pr	reventative maintenan	ice
Instrument	Known Standard	
Kit Used for Perfomance Check:	PrepFiler	PrepFiler BTA
Lot # Expiration	n Date	QC Date
Date Performance Check Completed		Analyst
Required records:	 QPCR res 	ults & setup worksheet
Results of Testing: Pass	Fail	(Choose One)
Comments:		
proval of COQC Record		
y Chief Criminalist, DNA am Leader, or designee:		Date*
		*If passed, this is the date released