



**Digest Buffer for DNA IQ**

**1. Required items:**

Tris-HCl, pH 7.5, 1M	1 ml
EDTA, 0.5 M	2 ml
NaCl, 5 M	1 ml
20% (w/v) SDS	2.5 ml
DI water	93.5 ml

**2. Instructions for Preparation:**

Mix the Tris-HCl, EDTA, NaCl, SDS, and DI water.

**3. Storage Conditions:**

Room Temperature

**4. Expiration Date:**

When the individual components expire, or twelve months after preparation, whichever comes first.

Lot # Tris-HCl 7.5/Exp. Date	Lot # EDTA/Exp. Date	Lot # NaCl/Exp. Date	Lot # SDS/Exp. Date	Amount Made	Date Made	Analyst	Solution Exp. Date