



**Leucomalachite Green (LMG)**

**1. Required items:**

Leucomalachite Green	0.05 g
Glacial Acetic Acid	33.5 ml
Zinc (mossy or powdered)	see instructions
DI water	16.5 ml

**2. Instructions for Preparation:**

Add Leucomalachite Green, Glacial Acetic Acid, and DI water, and dissolve.

**3. Storage Conditions:**

Refrigerate at 4°C in an amber bottle over Zinc

**4. Expiration Date:**

When the individual components expire, or twelve months after preparation, whichever comes first.

**5. QC needed.**

Lot# LMG / Exp. Date	Lot # Glacial Acetic Acid / Exp. Date	Lot # Zinc	Amount Made	Date Made	Analyst	Sol'n Exp. Date	QC Date